


Keysight Technologies

In Situ Investigations of Corrosion via SPM

Application Note

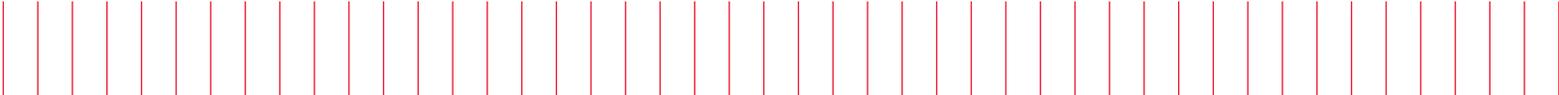

Introduction

Metals, minerals, and plastics are all subject to the destructive power of corrosion – affecting billions of dollars of repair and replacement costs per annum. It is not, however, a one-sided battle. Public and private sectors combat corrosion by pouring billions of dollars into research aimed at developing advanced corrosion-resistant materials, corrosion-resistant paints and coatings, and corrosion inhibitors.

Corrosion is an interfacial phenomenon, occurring at solid surfaces in contact with gases or liquids. Although it often has macroscopic consequences, corrosion typically begins at the atomic level. These factors make corrosion processes ideal candidates for study via *in situ* scanning probe microscopy (SPM), a highly sensitive surface technique that can be utilized in the presence of a wide range of corrosive solutions and gases. The versatility of *in situ* SPM enables direct, real-time observation of corrosional processes with atomic or near-atomic resolution, often providing valuable kinetic information as well as important structural data.

The use of *in situ* SPM as a key research tool for studying corrosion has been steadily gaining popularity over the past several years. Researchers have utilized SPM to investigate the structure of a variety of noble metal surfaces (1) and passivated surfaces (2, 3), electrochemical SPM to investigate the rate and mechanism of corrosion, and scanning tunneling microscopy (STM) to investigate the adsorption, orientation, and ordering of corrosion inhibitors on copper and silver surfaces (4). Information about the coverage and adherence of paints and other coatings is also easily accessible via use of these techniques.

Scanning probe microscopes, designed by Keysight Technologies, Inc., are uniquely suited to the *in situ* investigation of corrosion, providing a number of features designed specifically for these studies. For instance, Keysight's specially engineered electrochemical cell has a large volume-to-surface area ratio. The open design of the cell accommodates a large ring counter electrode, thus ensuring uniform distribution of current. Furthermore, removal of precipitates and gas bubbles can be accomplished by exchanging the solution in the cell without disrupting the experiment. Additionally, the Keysight environmental chamber can be used to control the amount of dissolved oxygen in the electrolyte or to introduce reactive gases. These features, combined with easy optical access, precision temperature control plates, and a truly modular design, make Keysight SPM systems an ideal choice for corrosion studies.

This application note describes three specific examples in which SPM has been used to study various aspects of metallic corrosion. More information about studies of corrosion using SPM can be found on Keysight's website or by contacting Keysight directly.

Corrosion of Steel

A study was conducted to investigate the effects of a sodium chloride solution on 304 stainless steel. In preparation, the sample was cut from a 304ss rod, polished to a roughness of less than 30 nm over 15 microns, and sonicated in deionized water for 20 minutes. The sample was transferred to the electrochemical cell and studied.

Figure 1(A) shows an image of the steel surface immediately upon immersion into 100 mM NaCl in purified, distilled water. The large inclusion in the bottom-left corner of the image was most likely a result of the polishing process. The sample was continually monitored while being held at open circuit in the corrosive solution. Images at 30 minutes (Figure 1(B)) and 150 minutes (Figure 1(C)) show the rapid dissolution of the inclusion and the corrosion of the surrounding stainless steel, suggesting that corrosion proceeds through a pitting mechanism.

Selective Dissolution of Aluminum

The tip-induced, selective dissolution of an aluminum foil is shown in Figure 2. The aluminum sample was imaged in 100 mM NaCl at high contact force, causing the formation of a recessed region. The dissolution mechanism is not simply rubbing away of the top layers of aluminum as one might assume from the image, as evidenced by the fact that the dissolution does not occur in the absence of chloride. Instead, the mechanism is thought to be a two-step process in which the passivating layer is first broken and then the chloride ions react with the underlying aluminum (5). Such experiments may represent a convenient method for studying the stability and strength of passivating coatings.

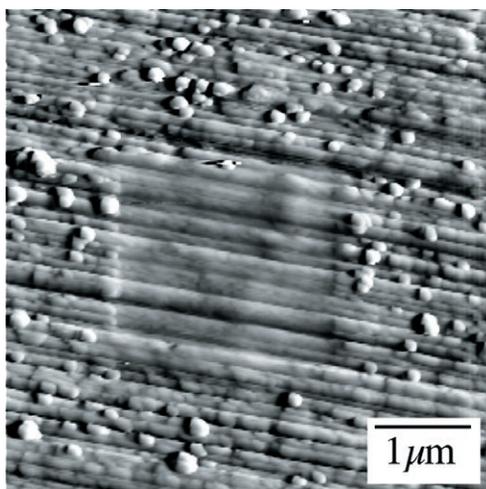


Figure 2. AFM image of the selective dissolution of aluminum in sodium chloride solution. The recessed region in the center of the image had been previously scanned at high force, accelerating the aluminum corrosion. Z range = 1 nm.

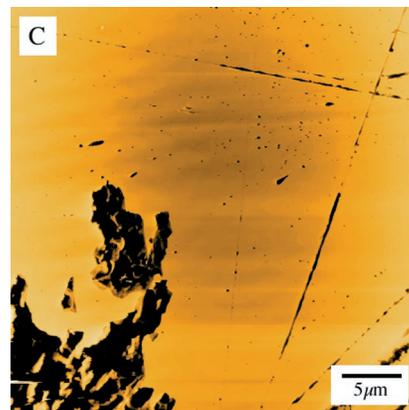
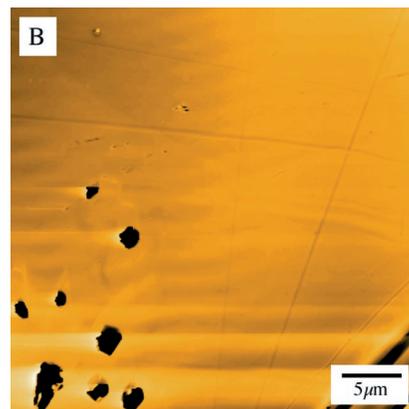
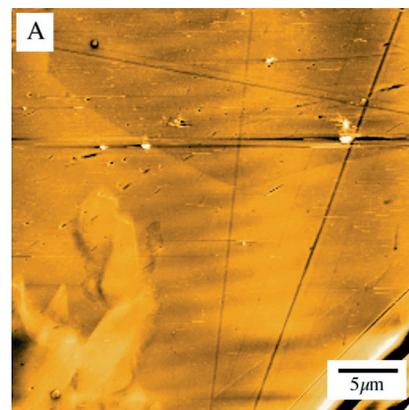


Figure 1. AFM images of the corrosion of 304 stainless steel in sodium chloride solution after (A) 4 minutes (z range = 50 nm), (B) 30 minutes (z range = 250 nm), and (C) 150 minutes (z range = 500 nm). Images courtesy of David Uhl, Arizona State University.

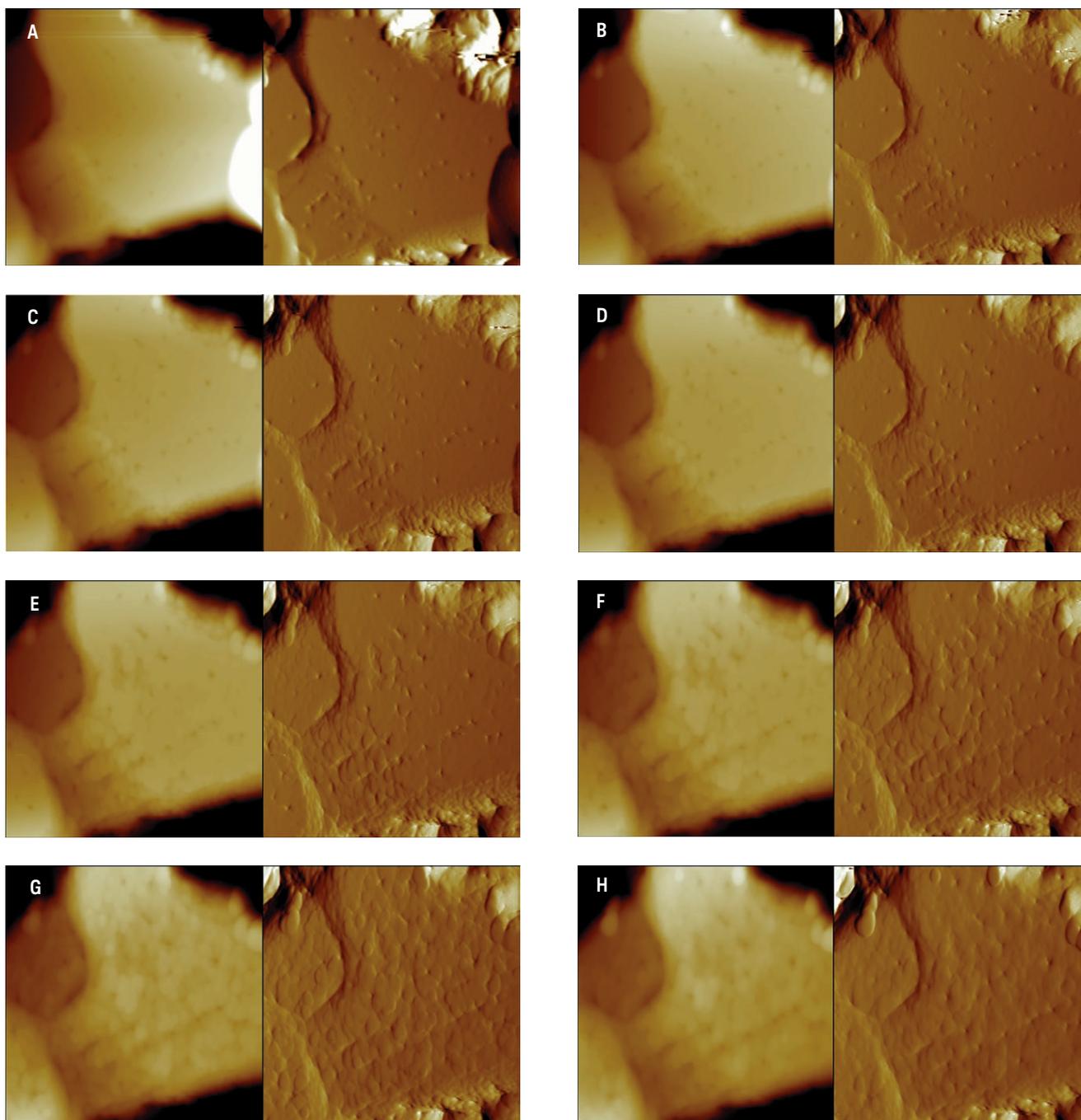


Figure 3. AFM height (left) and deflection (right) images of the corrosion of aluminum in 100 mM sulfuric acid after (A) 0 minutes, (B) 5.5 minutes, (C) 8 minutes, (D) 11 minutes, (E) 16 minutes, (F) 21.5 minutes, (G) 27 minutes, and (H) 43 minutes. Z range for height = 175 nm. Z range for deflection = 5 nm.

Kinetics of Aluminum Corrosion

Figure 3 shows a sequence of atomic force microscopy (AFM) images for the dissolution of an aluminum thin film in 100 mM sulfuric acid. In the preparation of the substrate, aluminum (99.999%, Puratronic) was deposited using physical vapor deposition onto mica in vacuum (base pressure = 5×10^{-9}). Imaging began immediately upon addition of 100 mM H_2SO_4 to the electrochemical cell; no efforts were made to eliminate oxygen from the electrolyte. The corrosion rate at open circuit was slow enough to capture a long image series that demonstrates the kinetics of aluminum corrosion, which proceeds through an apparent step dissolution process within terraces as well as at step sites and defects. Future studies will concentrate on the effects of corrosion inhibitors on the dissolution process.

References

1. Tao N.J., Pan J., Li Y., Oden P.I., DeRose J.A., Lindsay S.M., Surf. Sci., (1992), **272**, L338.
2. Maurice V., Yang W.P., Marcus P., J. Electrochem. Soc., (1994), **141**, 3016.
3. Maurice V., Talah H., Marcus P., Surf. Sci., (1993), **304**, 98.
4. Ge M., Zhong B., Klemperer W.G., Gewirth, A.A., J. Am. Chem. Soc., (1996), **118**, 5812.
5. Rouè L., Chen L., Guay D., Langmuir, (1996), **12**, 5818.

AFM Instrumentation from Keysight Technologies

Keysight Technologies offers high-precision, modular AFM solutions for research, industry, and education. Exceptional worldwide support is provided by experienced application scientists and technical service personnel. Keysight's leading-edge R&D laboratories are dedicated to the timely introduction and optimization of innovative and easy-to-use AFM technologies.

www.keysight.com/find/AFM

For more information on Keysight Technologies' products, applications or services, please contact your local Keysight office. The complete list is available at: www.keysight.com/find/contactus

Americas

| | |
|---------------|------------------|
| Canada | (877) 894 4414 |
| Brazil | 55 11 3351 7010 |
| Mexico | 001 800 254 2440 |
| United States | (800) 829 4444 |

Asia Pacific

| | |
|--------------------|----------------|
| Australia | 1 800 629 485 |
| China | 800 810 0189 |
| Hong Kong | 800 938 693 |
| India | 1 800 112 929 |
| Japan | 0120 (421) 345 |
| Korea | 080 769 0800 |
| Malaysia | 1 800 888 848 |
| Singapore | 1 800 375 8100 |
| Taiwan | 0800 047 866 |
| Other AP Countries | (65) 6375 8100 |

Europe & Middle East

| | |
|----------------|---------------|
| Austria | 0800 001122 |
| Belgium | 0800 58580 |
| Finland | 0800 523252 |
| France | 0805 980333 |
| Germany | 0800 6270999 |
| Ireland | 1800 832700 |
| Israel | 1 809 343051 |
| Italy | 800 599100 |
| Luxembourg | +32 800 58580 |
| Netherlands | 0800 0233200 |
| Russia | 8800 5009286 |
| Spain | 800 000154 |
| Sweden | 0200 882255 |
| Switzerland | 0800 805353 |
| | Opt. 1 (DE) |
| | Opt. 2 (FR) |
| | Opt. 3 (IT) |
| United Kingdom | 0800 0260637 |

For other unlisted countries:
www.keysight.com/find/contactus
 (BP-09-23-14)

This information is subject to change without notice.
 © Keysight Technologies, 2007 - 2014
 Published in USA, July 31, 2014
 5989-6895EN
www.keysight.com